



Glowing products for science

NucSpot® Nuclear Stains

Cell membrane-impermeant, nuclear-specific counterstains.
Suitable for fixed cells or staining dead cells in live cultures.



Product Description

NucSpot® Nuclear Stains are cell membrane-impermeant, nuclear-specific counterstains available in a variety of colors from green to near-infrared (near-IR).

- Nuclear-specific counterstains for fixed cells
- Selectively stains dead cells in live culture
- No wash required
- Color options for FITC, Cy@3, PE, Texas Red®, Cy@5, APC, Cy@5.5, and Cy@7 channels

Commonly used blue nuclear stains, such as DAPI and Hoechst, undergo photoconversion after exposure to UV wavelengths and cause cross-talk in other channels. NucSpot® Nuclear Stains were developed to address these photoconversion issues by offering bright and nuclear-specific counterstaining in a wide selection of colors from green to near-IR. They have minimal fluorescence until they bind to DNA and can be used for no-wash nuclear staining. Unlike other nucleic acid dyes such as propidium iodide (PI), TOTO®, TO-PRO®, and similar dyes that stain both the nucleus and cytoplasm, NucSpot® Nuclear Stains selectively stain the nucleus in fixed and permeabilized cells without the need for RNase treatment. NucSpot® Nuclear Stains also can be used for selective staining of dead cells in unfixed cell cultures for analysis by flow cytometry or fluorescence imaging. Several of the dyes can be continuously incubated with cells for multi-day imaging.

Call us : [800-304-5357](tel:800-304-5357)

Product attributes

Probe cellular localization	Nucleus
For live or fixed cells	For fixed cells
Assay type/options	Live/dead discrimination, Long term staining (24-72h), No-wash staining, Real-time imaging, Tissue staining
Detection method/readout	Fluorescence microscopy, Flow cytometry
Cell permeability	Membrane impermeant
Apoptosis/viability marker	Dead cell stain
Fixation options	Fix before staining (formaldehyde), Fix before staining (methanol), Permeabilize before staining
Colors	Green, Orange, Red, Far-red, Near-infrared
Storage Conditions	Store at 2 to 8 °C

NucSpot® Nuclear Stains

Product	Ex/Em	Detection Channel	Size	Catalog No.
NucSpot® 500/515 100 uL	497/513 nm 41040	FITC*	20 uL	41040-T
NucSpot® 555/570 100 uL	559/566 nm 41033	Cy@3 or PE*	20 uL	41033-T
NucSpot® 568/580 100 uL	572/583 nm 41036	Cy@3 or PE*	20 uL	41036-T
NucSpot® 594/615 100 uL	603/613 nm 41037	Texas Red® or PE-Texas Red®*	20 uL	41037-T
NucSpot® 650/665 100 uL	653/671 nm 41034	Cy@5 or APC*	20 uL	41034-T
NucSpot® 680/700 100 uL	683/707 nm 41035	Cy@5.5*	20 uL	41035-T
NucSpot® 750/780 100 uL	757/780 nm 41038	Cy@7 or APC-Cy@7*	20 uL	41038-T

* May show crosstalk in lower wavelength detection channels. Perform single-stain controls before combining with other probes.

† NucSpot® Far-Red is designed for flow cytometry optical systems and is not recommended for fluorescence microscopy.

Find the Right Dye for Your Application

Stain Dead Cells in Live Cultures

Exceptional Photostability

NucSpot® Nuclear Stains offer improved photostability over commonly used cyanine-based nuclear stains. The figure below shows how NucSpot® 650/665 is significantly more photostable than Thiazole Red Homodimer (TOTO® -3).

NucSpot® 500/515 is a green fluorescent nuclear stain for the FITC channel. In addition to bright nuclear counterstaining of fixed cells, it can also be used for multi-day live/dead staining of mammalian cells in culture. NucSpot® 500/515 is designed as an improved alternative to NucSpot® 470 which has limited use for cell viability tracking due to its instability in culture medium during long incubation periods.

NucSpot® 555/570 and NucSpot® 568/580 have orange and visible red fluorescence, respectively, and are nuclear-specific alternatives to PI and similar dyes.

NucSpot® 594/615 has deep red fluorescence for the Texas Red® channel.

NucSpot® 650/665 has far-red fluorescence with superior nuclear specificity compared to first-generation far-red nuclear stains such as RedDot™2 and Draq7™.

NucSpot® 680/700 and NucSpot® 750/780 are spectrally unique DNA stains for far-red and near-IR detection.

Biotium also offers [NucSpot® Far-Red](#), a flow cytometry stain developed as an improved alternative to 7-AAD. It shows less bleed-through fluorescence in the PE-Texas Red® channel compared to 7-AAD and is ideal for selective detection of dead cells or DNA content analysis by flow cytometry without RNase treatment.

Learn more about our full selection of [Cellular Stains](#) for the nucleus and other organelles.

TOTO and TO-PRO are registered trademarks of Thermo Fisher Scientific; Draq is a trademark of BioStatus Ltd; CY is a registered trademark of Cytiva.

This datasheet was generated on May 8, 2026 at 07:38:21 PM. Visit product page to check for updated information before use.

Product link: <https://biotium-woo.supremeclients.com/product/nucspot-nuclear-stains/>